



BioSurface Technologies Corporation

CDC Biofilm Reactor[®] (CBR) Operator's Manual



Prepared with help from the Center for Biofilm Engineering, Standardized Biofilm
Methods Laboratory for BioSurface Technologies Corporation

Contents

1. Purpose	3
2. CDC Biofilm Reactor® Description.....	3
3. CDC Biofilm Reactor® Standard Operating Procedure.....	4
A. Principle.....	5
B. Apparatus	5
C. Media and Reagents	9
D. General Instructions	9
a. Reactor Preparation.....	9
b. Batch Phase	10
c. Continuous Flow Phase	10
E. Sampling.....	10
a. Sampling Coupons	10
b. Disaggregation.....	11
c. Serial Dilution.....	11
d. Viable Colony Counts	11
F. Calculations	11
4. Repeatability	12
5. Troubleshooting.....	12
6. CDC Biofilm Reactor® Assembly	12
A. Stir Plate and Support Stand	12
B. Reactor Assembly	12
7. CDC Biofilm Reactor® Ancillary Equipment List	13
8. RTD Temperature Probe.....	14
9. RTD Temperature Probe Well Installation.....	15
10. Anaerobic CDC Biofilm Reactor® Assembly and Operation	16
11. CDC Biofilm Reactor® Reference Material	17
A. Standard Methods	17
a. ASTM (www.astm.org).....	17
b. USEPA (www.epa.gov).....	17
B. Training Videos.....	17
C. Publications	19

1. Purpose

The CDC Biofilm Reactor Operator's Manual is intended to serve as a guide for researchers interested in growing a laboratory biofilm under high shear conditions. It is the responsibility of the user to be familiar with basic microbiological concepts and techniques. Although a specific method is presented in the manual, the CDC Biofilm Reactor is suitable for modeling many different environments. Laboratory biofilms are engineered based upon the dynamics in the reactor where they grow. No one biofilm is better than any other, although one reactor may be a better choice for modeling a particular environment. It is the responsibility of the operator to choose which reactor best suits their research needs.

2. CDC Biofilm Reactor[®] Description

The CDC Biofilm Reactor (CBR, **Fig. 1**) is a one-liter vessel with an effluent spout at approximately 400 ml. Continuous mixing of the reactor's bulk fluid is provided by a baffled stir bar that is magnetically driven. An UHMW polyethylene top supports eight independent rods. Each rod houses three removable coupons (biofilm growth surfaces) for a total of 24 sampling opportunities. The coupons may be constructed from a variety of materials, for example: polycarbonate, porcelain, mild steel, stainless steel, PVC, vinyl, glass, etc. The coupons experience a consistent high shear from the rotation of the baffled stir bar. A researcher may vary shear by altering the baffle's rotational speed. The CBR operates as a continuous flow stirred tank reactor (CFSTR), meaning nutrients are continuously pumped into and flow out of the reactor at the same rate. A CFSTR offers the following advantages:

- the bulk fluid (all the fluid in the reactor) is assumed to be well mixed (no gradients)
- the surfaces within the reactor experience fluid shear (force opposing flow) from the mixing,
- the reactor can achieve steady state conditions which eliminates all terms that depend upon time in the mathematical representation of the reactor. This simplifies the mathematics necessary to calculate the reduction in biofilm accumulation as a result of a treatment,
- the residence time (time for one reactor sized volume of liquid to flow through the reactor) in the reactor is easily adjusted to select for biofilm growth for various species of bacteria,
- the nutrient feed may be specified and adapted for various bacteria or conditions (for example to represent a swimming pool environment or the white water associated with the paper industry),
- a portion of the reactor's effluent (waste) may be recycled back into the reactor or become the feed of a second reactor,
- with simple modifications, a CFSTR can be operated anaerobically (without oxygen). This increases the number of relevant environments that can be modeled using this design.

In addition to the list of advantages above, the CBR offers additional advantages:

- the biofilm growth surfaces (a.k.a. coupons) are interchangeable with other reactors that BioSurface Technologies manufactures and sells,
- 24 sampling opportunities,
- removable rods that allow for intermittent removal, control, and treated data from the same reactor, and
- a baffled stirrer provides for consistent shear across all the coupons.

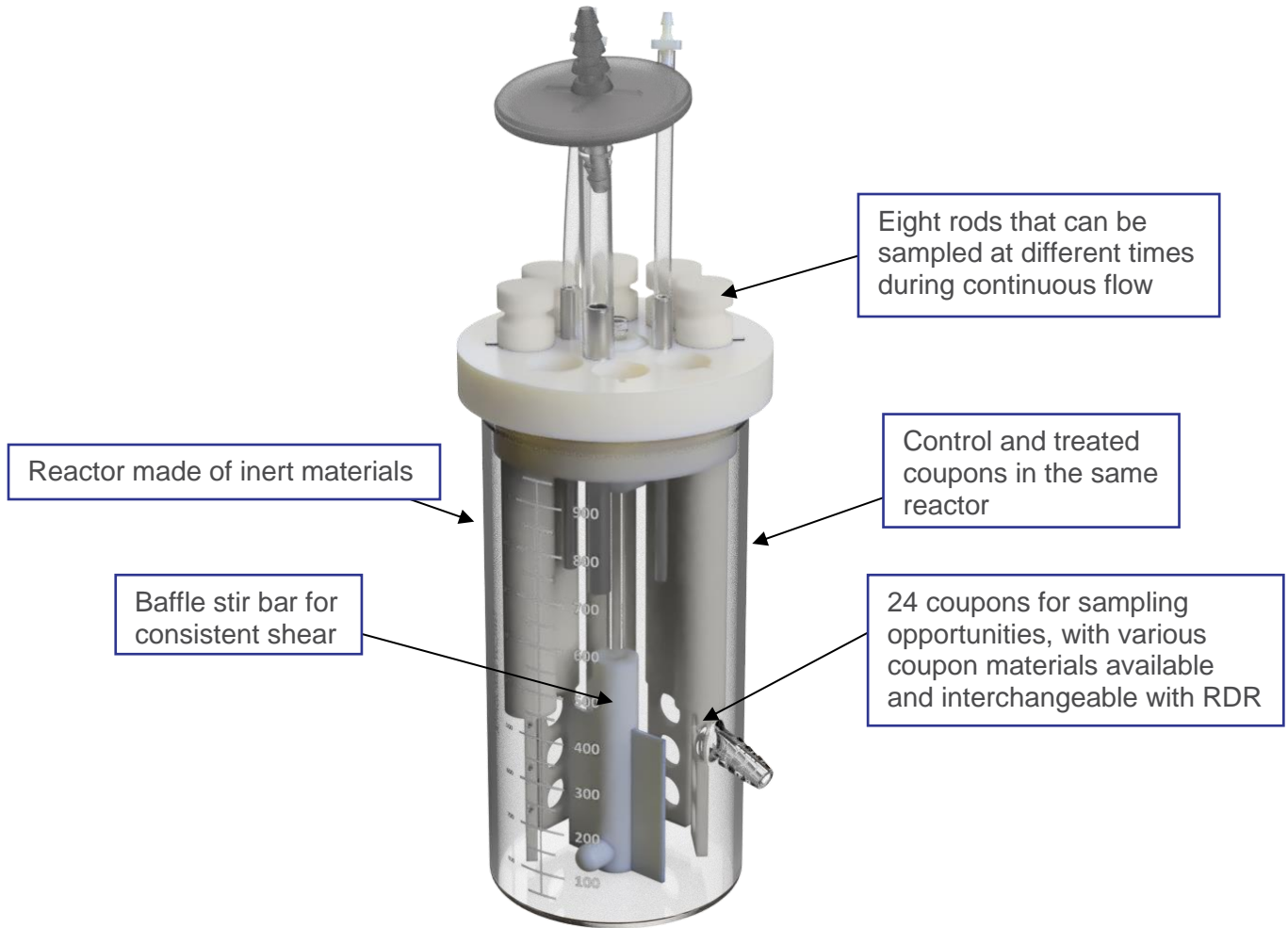


Figure 1: Picture of the CDC Biofilm Reactor.

3. CDC Biofilm Reactor® Standard Operating Procedure

The CBR was designed as a flexible reactor system, meaning that it can be easily adapted to model a variety of conditions in the laboratory. The following standard operating procedure (SOP) is just one suggested use. Operators are encouraged to modify the protocol to model the most appropriate biofilm for their research needs. A series of notes are included in the SOP that alert the operator to situations where special care must be taken.

Caution: All microorganisms should be handled according to safety recommendations for each individual species of microorganism. Decontamination of all media and equipment used during experimentation is required prior to disposal of media or re-use of equipment. It is the responsibility of the operator to inform themselves on these techniques.

A. Principle

Biofilm is defined as an accumulation of bacterial cells immobilized on a substratum and embedded in an organic polymer matrix of microbial origin. Biofilm bacteria exhibit a different phenotype from suspended bacterial cells of the same genotype. In this method, a laboratory biofilm is grown under high shear and continuous flow conditions. Biofilm accumulation is quantified by harvesting the biofilm from coupons of a known surface area, disaggregating the cells and polymer matrix, and performing viable plate counts.

B. Apparatus

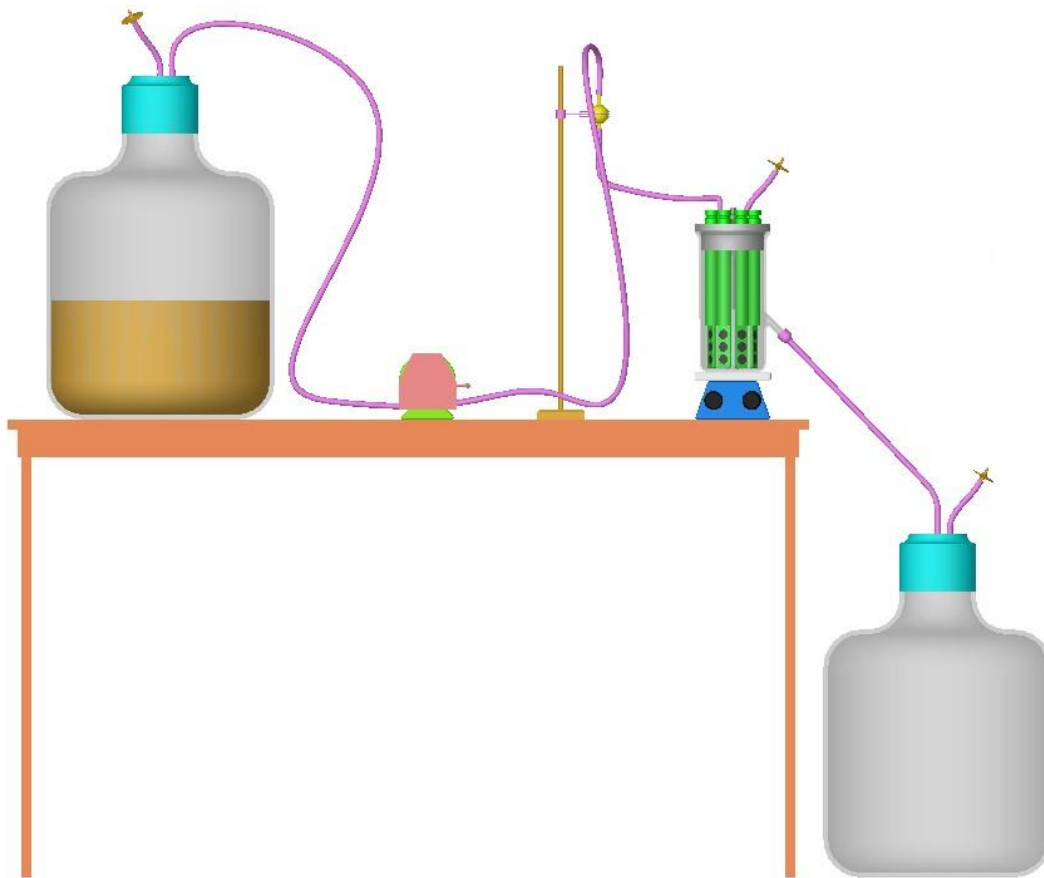


Figure 2a: Schematic of an assembled reactor system (Norris, 2003).



Figure 2b: Picture of an assembled reactor system (M. Mettler, CBE, 2020).

(a) Reactor Components

i) Borosilicate Glass Beaker: 1000mL beaker without conventional pour spout, 9.5 ± 0.5 cm diameter. Barbed outlet spout added so that bottom is at $400 \text{ mL} \pm 20\text{mL}$ mark. Angle the spout 30-45 degrees to ensure drainage. Spout should accommodate silicon tubing with ID of 8-11 mm.

NOTE 1: The rods (B.a.iii) and baffle (B.a.v) will displace approximately 50 mL of liquid when the system is completely assembled. Therefore, an outlet spout at the 400 mL mark will result in a volume of approximately 350 mL during operation. Before beginning an experiment, confirm the actual reactor operating volume when the reactor is fully assembled.

ii) Reactor top, Figure 2: UHMW polyethylene top equipped with 3 holes accommodating 6-8 cm long pieces of stainless steel or other rigid autoclavable tubing with OD of 5-8 mm for media inlet, air exchange and inoculation port. Center hole, to accommodate the glass tubing used to support the baffle assembly (B.a.v). Eight rod holes notched to accommodate stainless steel rod alignment spike. Rubber gasket, to fit between reactor top and reactor beaker.

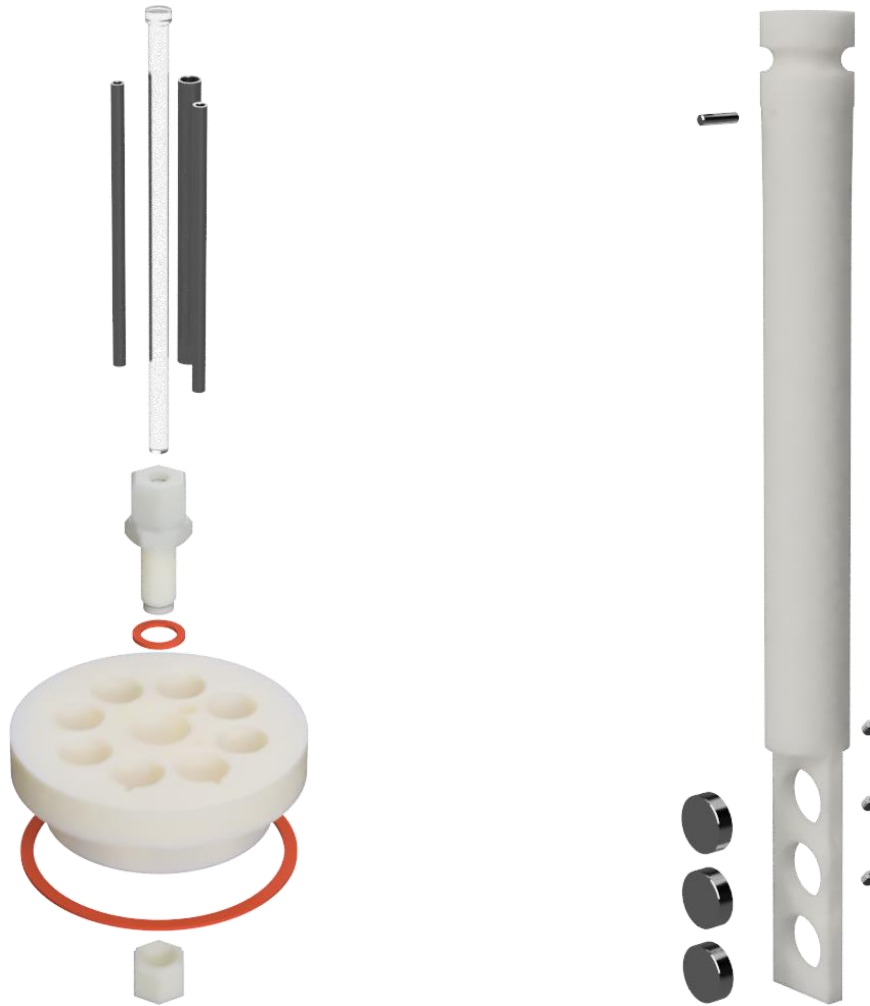


Figure 3a (Left): Expanded schematic of reactor top.
Figure 3b (Right): Expanded schematic of coupon rod and coupons.

iii) Polypropylene rods, Fig. 3b: Eight polypropylene rods, machined to hold three coupons (B.a.iv) at the immersed end. 316 Stainless Steel set screws imbedded inside to hold coupons in place. Rods fit into holes in reactor top and lock into preformed notches.

iv) Polycarbonate coupons, Fig. 3b: Provide a surface for biofilm growth. Twenty-four cylindrical polycarbonate coupons with a diameter of 1.27 cm, thickness of approximately 3.0 mm.

v) Stir Blade Assembly (Baffle), Fig. 4: To create constant shear/mixing. PTFE blade fitted into cylindrical PTFE holder and held in place with a magnetic stir bar. PTFE holder fits onto a glass rod fitted into the reactor top (Fig. 3a). The glass rod is held in place with a compression bulk-head fitting and acts as a support for the moving blade assembly.

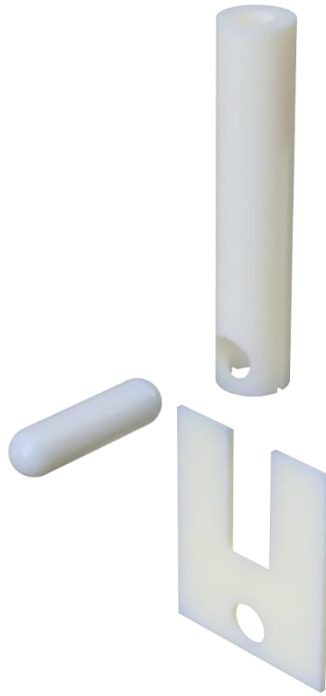


Figure 4: Expanded schematic of baffled stir bar.

- (b) *Glass Flow break*: Autoclavable and accommodating tubing of ID 3.2 mm. Insert vertically into media inlet tube to prevent contamination from reactor to media carboy.
- (c) *Silicone tubing*: Two sizes of tubing, one with ID 3.2 mm (port OD=4.75 mm) and the other with ID 7.9 mm (port OD=7.94 mm). Both must withstand sterilization. Use to connect media carboy to reactor inlet; beaker outlet to waste carboy; air vent to reactor top; media carboy air vent.
- (d) *Carboy*: Two 20L autoclavable carboys for media and waste.
- (e) *Carboy Tops*: Two threaded tops each equipped with 2 barbed fittings to accommodate tubing ID 3.1 mm for nutrient or waste line and air vent attachment.
- (f) *Magnetic Stir plate*: Capable of operation at 100-400 rpm. A digital stir plate is recommended.
- (g) *Pipette*: Capable of accurately dispensing 1.0 mL.
- (h) *Micropipette*: Capable of accurately dispensing 0.01 mL.
- (i) *Top Loading Balance*: For weighing reagent ingredients; sensitive to 0.01 gm.
- (j) *Homogenizer Probe*: Capable of mixing at $20,500 \pm 5000$ rpm in a 5-10 mL volume and able to withstand autoclaving or other means of sterilization.
- (k) *Vortex*: Any vortex mixer that will ensure proper mixing of culture tubes.
- (l) *Peristaltic Pump*: For pumping media into reactor during continuous flow phase; pump head capable of holding selected tubing (accommodates tubing ID/OD).
- (m) *Wooden Applicator Sticks*: Used for scraping biofilm from coupon surface. Sterile.
- (n) *Hemostat*: Stainless steel hemostat clamp with curved tip to remove and hold coupon in place while scraping. Use after flame sterilizing.
- (o) *Inoculating Loop*. Used to introduce bacteria into inoculum flask. Sterile.
- (p) *Bunsen or Alcohol burner*: For flame sterilizing inoculating loop and hemostat.
- (q) *Culture Tubes with caps*: Any with a volume capability of 10 mL and diameter no less than 6 cm. Recommended size is 16 x 125mm borosilicate glass with threaded opening.
- (r) *Sterilizers*: Any steam sterilizer capable of producing the conditions of sterilization.
- (s) *Colony Counter*: To aid in counting plates.
- (t) *Clamp Stand and Clamp*: To hold glass flow break and media inlet tube vertically above reactor top. Clamp with 0.5 cm minimum grip size.
- (u) *Petri Dish*: 100 x 15mm, plastic, sterile, for transporting rods from reactor to workstation.

(v) *Plexiglas Board*: Or other surface, used as a sampling platform on which to hold and scrape samples. Use after disinfecting with 70% EtOH.

(w) *Environmental Shaker*: Capable of maintaining temperature of 37°C.

C. Media and Reagents

NOTE 2: All reference to water as diluent or reagent means distilled water or water of equal purity.

(a) *Inoculum culture media*: 100mL of 300mg/L soybean-casein digest media or an equivalent general bacterial growth medium. Tryptic soy broth is recommended. Sterilize for 20 minutes. Inoculate media with a single colony from a bacterial isolation plate using sterile inoculating loop. *Pseudomonas aeruginosa* (ATCC 700888) is the bacteria used in this method. Incubate at $35 \pm 2^\circ\text{C}$ while shaking for 18-24 h.

(b) *Batch culture media*: 500mL of 300mg/L TSB. Sterilize in reactor for 20 minutes.

(c) *Continuous flow media*: 20L of 100mg/L TSB. Sterilize 20L of water then aseptically add concentrated broth (2.0 gm TSB/500mL, sterilized 20 minutes) so that the final volume equals 20 L and final concentration equals 100 mg/L.

(d) *Buffered Dilution Water*: 0.0425 g/L KH_2PO_4 distilled water, filter sterilized and 0.405 g/L $\text{MgCl}\cdot 6\text{H}_2\text{O}$ distilled water, filter sterilized. Aseptically fill sterile, capped test tubes with 9 mL of sterile dilution water.

(e) *Bacterial Plating Medium*: R2A agar is recommended.

(f) *95% Ethanol*: For dipping hemostat prior to flame sterilizing.

(g) *70% Ethanol*: For rinsing homogenizer probe and general bench top clean up.

D. General Instructions

a. Reactor Preparation

i) Clean reactor coupons by immersing in a 1:1000 dilution of laboratory soap and sonicating for 3-5 minutes. Rinse with tap water until no more soap appears then sonicate again in reagent grade water. A final 2-hour soak in 2M HCl is recommended then a final rinse with reagent water (**do not soak hydroxyapatite or other acid susceptible coupons in HCl; NOTE 3**). Let coupons air dry. Place a coupon into each hole of the reactor rod, leaving the top of the coupon flush with the inside rod surface (sampled surface). Hold in place by tightening the set screw adjacent to the hole.

NOTE 3: Check for material compatibility before soaking coupons in 2 M HCl. Some coupon materials, such as hydroxyapatite, carbon steel, or stainless steel, should not be soaked in 2M HCl. If the material is incompatible with this acid, investigate alternative cleaning solutions.

NOTE 4: Visually inspect coupon surface topography for any deformations before using. Discard damaged coupons.

ii) Assemble all reactor parts including rods, coupons, baffle, bacterial air vent and tubing. Air vent and tubing are attached to the rigid tubes in the reactor top (B.a.ii). Use a small piece of flexible tubing to attach the air vent to the reactor top. Check assembly on stir plate to ensure baffle rotates without scraping rods. Clamp effluent tube. Foil the exposed ends of all tubing.

iii) Add 500 mL non-sterile batch culture media (C.b). Sterilize reactor system using an autoclave for 20 minutes at 121°C.

NOTE 5: Autoclaving longer than 20 minutes could cause unnecessary degradation to the reactor or its components and is not advised. The reactor and its components have not been approved for temperatures above 121°C, and any damage caused by a higher temperature will not be covered under warranty.

b. Batch Phase

- i)* Place reactor on stir plate. Be sure gases may pass freely through the bacterial air vent and effluent tubing is clamped shut.
- ii)* Aseptically inoculate the reactor by injecting 1mL of previously prepared inoculum culture (C.a) into the batch media through an inoculum port in reactor top.

NOTE 6: The viable bacterial density of the inoculum should equal 10^8 cfu/mL. Confirm this number by diluting and plating a sample from the inoculum flask.

- iii)* Start the baffle rotating at 125 ± 5 rpm.

NOTE 7: The speed at which the baffle rotates directly determines the amount of shear stress that the biofilm experiences. Ruggedness testing showed that biofilm accumulation on the coupons is sensitive to changes in the baffle's rotational speed. ***The baffle rotational speed is a critical factor that must be controlled. If a digital stir plate is not available, then use a strobe light to confirm rotational speed.***

- iv)* Operate the reactor in batch phase for 24 hours.

c. Continuous Flow Phase

- i)* Aseptically connect the nutrient tubing line to the carboy containing the continuous flow nutrient broth (C.c).
- ii)* Remove foil and place the end of the effluent tubing into a waste carboy, unclamp.
- iii)* Pump a continuous flow of media into the reactor at a flow rate equal 11.67 ± 0.2 ml/min.

NOTE 8: Flow rate is calculated by dividing the reactor volume by the residence time. The residence time is 30 minutes. The reactor volume is approximately 350 ml (**NOTE 1**). The operator should set an exact flow rate based upon the measured fluid volume in the reactor when the baffle, rods, and coupons are in place to achieve an exact 30-minute residence time. The reactor residence time is a critical parameter that is specific to the bacterial species used during the experiment. To select for biofilm growth in the reactor, the residence time must be less than the doubling time for the suspended cells. This will result in the suspended cells washing out of the reactor, leaving only biofilm.

- iv)* The reactor is operated with continuous flow for 24 hours.

E. Sampling

a. Sampling Coupons

- i.)* Set-up sampling materials: vortex, homogenizer, culture tubes, pipettes, wooden sticks, flame sterilized hemostat, sampling platform.
- ii.)* Turn off stir plate and pump.
- iii.)* Remove a single rod by pulling straight up. Collect any drips in a sterile petri plate held beneath rod.

- iv.) Loosen set screw to release coupon and remove using flame sterilized hemostat, being careful not to disturb the biofilm on coupon surface that will be scraped.
- v.) Hold coupon in place on a disinfected Plexiglas Board with hemostat and scrape the side of the coupon that faced the baffle with a sterile wooden stick. Hold the stick perpendicular to the surface when scraping.
- vi.) Remove cap of first 9 mL dilution tube and rinse stick by swirling and tapping on bottom of tube. Repeat scraping and rinsing steps 3 to 4 times.
- vii.) Hold coupon above dilution tube and rinse scraped surface with 1 mL dilution water. The total volume in the tube equals 10 mL. Record this number as the volume scraped into **Eq. 1**.

b. Disaggregation

- i.) Homogenize the dilution tube containing sample at 20,500 rpm for 30 s using sterile homogenizer probe.
- ii.) Clean the probe between samples by rinsing for 30 s at 20,500 rpm in a sterile dilution blank followed by a rinse at 20,500 rpm for 15 s in 75% ethanol. Let the probe soak in ethanol for 1 min. Rinse probe two more times with dilution blanks at 20,500 rpm for 30 s each.

NOTE 9: Homogenization breaks up biofilm clumps to form a homogeneous cell suspension. Improper disaggregation will result in an underestimation of viable cells present in the sample.

c. Serial Dilution

- i.) Serially dilute the disaggregated biofilm sample according to standard microbiological techniques.

d. Viable Colony Counts

- i.) Plate each dilution in duplicate on R2A nutrient agar. Any standard plating technique, such as spread plating or drop plating, is acceptable.
- ii.) Incubate plates at 35 ± 2°C for 18-24 hours.
- iii.) Count the appropriate number of colonies for the plating method used.

F. Calculations

a) Calculate the log density for one coupon using **Equation 1**:

$$LOG_{10}\left(\frac{cfu}{cm^2}\right) = LOG_{10}\left[\left(\frac{\left(\frac{mean\ cfu}{plate}\right)}{vol.\ of\ sample\ plated}\right) \times \left(\frac{volumescraped\ in\ to}{surface\ area\ scraped}\right) \times (dilution)\right]$$

b) The mean biofilm density is the average of the individual coupon log densities.

4. Repeatability

For the CBR SOP, the estimated within-experiment variance is 0.1980 and the estimated between-experiment variance is 0.1461. The estimated repeatability standard deviation is 0.5866 of which 58% is attributable to within-experiment variation and 42% is attributable to between-experiment variation. This repeatability standard deviation pertains to a protocol that samples only one coupon per experiment. The repeatability standard deviation for a protocol that requires sampling n coupons per experiment is $\sqrt{[(0.1980 / n) + 0.1461]}$. For example, if the protocol specifies $n=3$ coupons and the reported log density is the mean of the three individual coupon log densities, then the repeatability standard deviation is 0.46 of which 31% is attributable to within-experiment variation and 69% is attributable to between-experiment variation.

5. Troubleshooting

1. Always check the alignment of the rubber gasket (**Fig. 3a**).
2. Visually confirm that the inoculum reaches the reactor's bulk fluid.
3. The top of the glass rod is the end with the expanded head (**Fig. 3a**). This design holds the baffle in the center of the reactor while allowing it to rest on the bottom for proper mixing.
4. Tighten coupon set screws before the reactor is autoclaved to ensure correct alignment.
5. If a rod pops out of the reactor during autoclaving, immediately push it back into the reactor top and continue with experiment. If the operator suspects the reactor's sterility has been compromised, start over from the beginning of the protocol.
6. Always ensure that the reactor's effluent spout allows for proper draining.
7. If rod alignment peg falls out (**Fig. 3b**), replace before beginning an experiment.
8. Confirm that the bacterial air vent is working properly before an experiment begins, so that the reactor can properly exhaust during autoclaving.

6. CDC Biofilm Reactor[®] Assembly

NOTE 10: The CDC Biofilm Reactor System (CBR 90) comes with and without a stir plate and support stand. These assembly directions assume a stir plate and support stand are used.

A. Stir Plate and Support Stand

- i. Fix the aluminum rod into the back of the stir plate in the support clamp.
- ii. Mount the support ring in a clamp and attach to the support rod. Mount the ring low on the stand near the top of the stir plate.
- iii. Mount the 3-prong clamp onto the support rod with a clamp. Mount the 3-prong clamp high on the support rod.

B. Reactor Assembly

- i. Nutrient Supply Tubing: Attach tubing from your media reservoir to the glass flow break inlets.
- ii. Attach effluent tubing. Use minimum 1/4"(6 mm) ID and direct the effluent tubing to the drain or collection reservoir.
- iii. Insert the coupons into the coupon rods, flush with the inner rod surface.
- iv. Insert the rotor into the glass reactor vessel.
- v. Mount the reactor lid onto the glass vessel while inserting the glass rod into the center of the blade rotor.
- vi. Insert the glass vessel bottom through the stand ring support. Adjust the location of the ring support to center the glass vessel in the middle of the stir plate.

- vii. Adjust the 3-prong support clamp to hold the influent tubing and glass flow breaks above the reactor vessel.
- viii. Insert rod coupon holders into the reactor vessel through the top access ports. Turn the coupon holders so the face of the coupons is perpendicular to the blade on the vane rotor (stainless steel alignment pin aligns with slot on the reactor lid) and push the coupon holders down into the ports.

7. CDC Biofilm Reactor® Ancillary Equipment List

NOTE 11: *BioSurface Technologies does not sell or supply the ancillary equipment described below. This or comparable equipment is required to complete the reactor system set-up. Please check with your local vendors for availability and current pricing. Equipment listed is not recommended equipment, but an aid to help you identify compatible equipment.*

Pump: There are many types of peristaltic pumps available, and you may opt for alternative types and pump head configurations (multi-channel versus single channel, higher/lower rpm range, etc.). Depending on what you need to add to the reactor during operation, you may require more than 1 pump or a multi-head pump (media for growth and biofilm treatment chemical for some duration that may require an additional pump).

- **Masterflex® L/S® Digital Precision Modular Drive with Remote I/O and Benchtop Controller, 1 to 100 rpm; 90 to 260 VAC**
(VWR Item#: MFLX07557-10)
- **Masterflex® L/S® Easy-Load® II Pump Head for Precision Tubing, PPS Housing, SS Rotor**
(VWR Item#: MFLX77200-60)

Tubing: The tubing you choose depends on the chemical compatibility, gas permeability, wear resistance in peristaltic pumps, and pricing. You must choose the tubing that best fits your needs. C-Flex tubing (listed below) is similar to silicone tubing but has a low gas-permeability compared to silicone. If gas-permeability is not an issue, standard silicone tubing is acceptable.

- **Masterflex® L/S® Precision Pump Tubing, C-Flex®, L/S 16; 25 ft**
(VWR Item#: MFLX06424-16)
 - Will fit the 3/16" barbed port fittings on the top of the CBR and connect to the media supply reservoir.
- **Masterflex® L/S® Precision Pump Tubing, C-Flex®, L/S 18; 25 ft**
(VWR Item#: MFLX06424-18)
 - Used to connect the effluent port (3/8" barbed port) on the reactor to the spent media collection vessel.
- **Masterflex® Transfer Tubing, C-Flex®, Opaque White, 1/4" ID x 7/16" OD; 25 Ft**
(VWR Item#: MFLX06424-72)
 - A few lengths and adapters to get from the carboy to the smaller diameter tubing, and as a siphon tube inside the carboy.
- **Masterflex® Fitting, Nylon, Straight, Hose Barb Reducer, 1/4" ID x 1/8" ID; 10/PK**
(VWR Item#MFLX30622-28)
 - Needed to get from the 3/16" or 1/4" ID to the 1/8" tubing.

Carboy: Carboys should be selected based on experiment needs and may be larger or smaller than what is suggested below. Ported lids can be purchased from suppliers, but standard lids are easily converted to ported lids using the following fittings or similar.

- **Azlon® Bottle, Rounded Octagonal, Polypropylene, Dynalon, 10L**
(VWR Item#: 30620-188)
- **Azlon® Bottle, Rounded Octagonal, Polypropylene, Dynalon, 20L**
(VWR Item#: 76210-720)
- **Nalgene® Barbed Bulkhead Fittings, Thermo Scientific, 6.4 mm (1/4")**
(VWR Item#: 16331-102)
- **Nalgene® Barbed Bulkhead Fittings, Thermo Scientific, 12.7 mm (1/2")**
(VWR Item#: 16225-232)
- **Cole-Parmer PTFE Syringe Filters, Non-Sterile; 0.45 µm, 50 mm Diameter**
(Cole Parmer P/N: EW-02915-30)

Suggested Ancillary Equipment Suppliers:

VWR: 800-932-5000 (www.vwr.com)

Cole Parmer: 800-323-4340 (www.coleparmer.com)

Fisher Scientific: 800-766-7000 (www.fishersci.com)

8. RTD Temperature Probe

(for use with the digital stir/hot plate)

The CDC Biofilm Reactor with Stir/Hot Plate is provided with an RTD Temperature Sensor for temperature control of the Stir/Hot Plate. The standard part provided for the RTD Sensor is an RTD sensor well (P/N CBR 2218). The location of the temperature sensor in the reactor vessel is critical to providing an accurate temperature measurement of the liquid media within the reactor. This well consists of nylon, 316 stainless steel, or anodized aluminum tubing sealed at the bottom. The well allows insertion of the sensor into the reactor vessel but does not directly contact the liquid. This location of the RTD Sensor allows accurate temperature control but reduced temperature change response times. Allow extra time for the stir/hot plate to bring the temperature within the vessel to the desired set-point. Remove the sensor from the well prior to autoclaving the vessel.

NOTE 12: To increase accuracy of the temperature readings, the well should be partially filled with mineral oil to allow for better temperature conductivity to the probe inside the well.

ALTERNATE RTD SENSOR LOCATIONS

T-Fitting on Effluent Port: The RTD sensor tip can be placed in contact with the effluent flow by using a t-fitting to attach to the effluent port of the glass vessel. At higher temperature operations, the sensor may need to be inserted into the reactor vessel to provide accurate temperature control. The RTD sensor may be inserted into the effluent port using the t-fitting (inserted the length of the base of the tee, with the effluent flow exiting to the waste collection vessel through the perpendicular port on the t-fitting; make sure the fitting, sensor seal, and vessel port are liquid tight prior to operation).

To create this t-fitting setup for effluent flow temperature measurement, use a 1/2-inch barbed t-fitting (preferably nylon or another autoclavable plastic). Cut a short section of 1/2-inch silicone tubing and attach to one side and the top of the fitting. Cut another length of 3/8-inch OD silicone tubing and insert into the top of the 1/2-inch tubing until it bottoms out. Cut another length of 1/4-inch silicone tubing and slide it onto the temperature probe, and then insert the

probe with tubing into the top of the t-fitting. This stacked-tubing method should create a good seal to prevent leaking. The entire fitting can then be attached to the effluent barbed port.

Through a Coupon Holder Rod Port: The RTD sensor may be mounted through one of the top ports (coupon holder rod ports) using a properly sized tapered hole stopper, bored to fit the temperature probe. This location allows direct contact of the reactor media to the temperature sensor and will provide the most accurate temperature control.

9. RTD Temperature Probe Well Installation (for use with the digital stir/hot plate)

The central glass shaft can be replaced with an RTD sensor well (P/N CBR 2218). This well consists of rigid nylon, 316 stainless steel, or anodized aluminum tubing sealed at the bottom. The well allows insertion of the sensor into the reactor vessel but does not directly contact the liquid. This location of the RTD sensor allows accurate temperature control, but reduced temperature change response times.

To replace the glass shaft with the temperature well, two 1 1/16" sockets (6-point, 18 mm) and socket drive are required. Place one of the sockets in a vise, pointing up. Insert the CDC lid with glass shaft into the immobilized socket and align the hex nut on the bottom side of the lid into the hex socket. Place the other socket onto the hex nut on the top of the lid. Using the socket driver, turn the nut and lid counterclockwise to loosen the bottom nut. Remove the bottom nut (if the top nut loosens before the bottom nut, place a flat bladed screw driver into the lid recess next to the compression fitting holding the glass rod and apply pressure to the screw driver to prevent the compression fitting from rotating, while turning the lid counter-clockwise. This rotation will loosen the compression nut on the bottom of the lid.). Press the bulkhead compression fitting out of the lid (the bulkhead compression fitting is not threaded to the lid). Insert the new bulkhead fitting with sensor well and compression gasket installed into the lid and install the bottom hex nut. Place the lid onto the socket held in the vise and tighten the top hex nut using the socket and driver just until the top bulkhead gasket is compressed (remove the top socket, look down into the compression fitting recess. The compression gasket will start to bulge out when it is compressed. Over-compression will cause the gasket to bulge on one side significantly and un-seat the gasket from between the two mating surfaces).

The RTD sensor well should now be ready for use. The RTD sensor will slide into the well from the top of the reactor. The RTD sensor does not directly contact the liquid in the vessel, which reduces its response time. Allow extra time for the stir/hot plate to bring the temperature within the vessel to the desired set-point. Remove the sensor from the well prior to autoclaving the vessel.

10. Anaerobic CDC Biofilm Reactor[®] Assembly and Operation

Do Not Pressurize or Operate Under Vacuum

The CDC Biofilm Reactor lid modification for anaerobic gas containment does not make the CDC Biofilm Reactor a pressure or vacuum containment vessel. The seals added to the lid are to aid in containment of anaerobic gases and to reduce oxygen migration into the vessel during anaerobic microbial growth. The CDC Biofilm Reactor vessel should always be vented.

The rubber flange gasket on the underside of the reactor lid can be torn easily. Ensure that the setscrews on the coupon holder rods are not protruding from the sides of the holder rods before inserting them into the reactor lid.

BioSurface Technologies does not recommend autoclaving with the lid clamps in place, nor is this necessary. All components of the Anaerobic CDC Biofilm Reactor (CBR 90-3) are autoclavable and re-useable.

Assembly Instructions

The modified lid system includes an O-ring seal and compression flanges, and a silicone flange gasket on the underside of the reactor lid. The silicone flange gasket will provide a light seal around each coupon holder rod when they are fully inserted into the lid. The compression rings and silicone O-ring, along with four stainless steel wingnuts, create a seal between the lid and the glass vessel.

The bottom compression ring (with screws mounted in four holes) should be lowered over the top of the glass vessel so that the screws are facing up. The plastic split ring is installed over the top of the glass vessel by opening the ring slit and working the ring over the glass vessel lip. The silicone O-ring is installed around the top of the glass vessel near the vessel lip. The top compression ring is placed over the top of the CDC Biofilm Reactor lid after it is installed, and the 4 holes are aligned with the screws on the bottom compression ring. The 4 wingnuts are then used to compress the rubber O-ring and underside of the reactor lid (avoid over-tightening the wingnuts by tightening them evenly around the lid to a finger-tight fit).

The lid ports on the Anaerobic CDC Biofilm Reactor have been replaced with polymer tubing. The coupon holder rod setscrews are 316 stainless steel and are not compatible with all anaerobic gases or anaerobic conditions (they will corrode under H₂S or acidic anaerobic conditions). A set of nylon setscrews has been provided as replacement items for conditions which are not compatible with stainless steel.

A clamp or containment ring should be used when mounting the reactor on a stir plate to prevent the reactor from tipping (the anaerobic containment system makes the reactor somewhat top-heavy).

11. CDC Biofilm Reactor® Reference Material

A. Standard Methods

a. ASTM (www.astm.org)

ASTM E2562

Standard Test Method for Quantification of *Pseudomonas aeruginosa* Biofilm Grown with High Shear and Continuous Flow using CDC Biofilm Reactor

ASTM E2871

Standard Test Method for Evaluating Disinfectant Efficacy against *Pseudomonas aeruginosa* Biofilm Grown in CDC Biofilm Reactor using Single Tube Method

ASTM E3161

Standard Practice for Preparing a *Pseudomonas aeruginosa* or *Staphylococcus aureus* Biofilm using the CDC Biofilm Reactor

b. USEPA (www.epa.gov)

MB-19

Preparing a *Pseudomonas aeruginosa* or *Staphylococcus aureus* Biofilm using the CDC Biofilm Reactor.

MB-20

Single Tube Method for Determining the Efficacy of Disinfectants against Bacterial Biofilm

B. Training Videos

Center for Biofilm Engineering Standardized Biofilm Methods Laboratory. "Single Tube Method Video." *YouTube*, uploaded by BiofilmFirst, 20 Aug. 2018, www.youtube.com/watch?v=vwpkxOqYkhU.

---. "Intro for Standardized Biofilm Methods Laboratory Training Video." *YouTube*, uploaded by BiofilmFirst, 10 Oct. 2016, www.youtube.com/watch?v=Rxi6pn--JOA.

---. "Chapter 01 Cleaning reactor components." *YouTube*, uploaded by BiofilmFirst, 26 July 2016, www.youtube.com/watch?v=3eCeOyDASzk.

---. "Chapter 02 Cleaning Coupons." *YouTube*, uploaded by BiofilmFirst, 26 July 2016, www.youtube.com/watch?v=7AF21JBxB9w.

---. "Chapter 03 Baffle Assembly." *YouTube*, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=vbp50MsEj_Y.

---. "Chapter 04 Bacterial Air Vent Assembly." *YouTube*, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=SVHvfEbH6e8.

---. "Chapter 05 Inoculation Port Assembly." *YouTube*, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=GcYPbg3J42U.

---. "Chapter 06 Continuous Flow Tubing Assembly." *YouTube*, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=mbLzCe0DaDo.

---. "Chapter 07 Waste Carboy Assembly." YouTube, uploaded by BiofilmFirst, 25 July 2017, www.youtube.com/watch?v=qiTrgdJ9DPg.

---. "Chapter 08 Coupon Installation." YouTube, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=Aclmg_H2xvE.

---. "Chapter 09 Measuring Reactor Volume." YouTube, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=-RXF2JBhZWg&t=3s.

---. "Chapter 10 Calculating Pump Flow Rate." YouTube, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=igMfU8P04vQ.

---. "Chapter 11 Continuous Flow Media Carboy Assembly." YouTube, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=LnKcWwEhSgU.

---. "Chapter 12 Waste Carboy Assembly." YouTube, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=yWiNLZZR-4o.

---. "Chapter 13 Preparing Reactor and Medium for Sterilization." YouTube, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=zIIYha4_c-s.

---. "Chapter 14 Addition of Batch Medium to the Reactor." YouTube, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=xtshqA66BU8&t=1s.

---. "Chapter 15 Inoculation of the Reactor and Start of Batch Phase." YouTube, uploaded by BiofilmFirst, 17 Aug. 2016, www.youtube.com/watch?v=GHUf4Em4JEs.

---. "Chapter 16 Preparation of Continuous Flow Medium." YouTube, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=YsiiyWSd9ZI.

---. "Chapter 17 End of Batch Phase and Start of Continuous Flow Phase." YouTube, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=NfHZ0BwGAuY.

---. "Chapter 18 End of Continuous Flow Phase and Sampling Day." YouTube, uploaded by BiofilmFirst, 27 July 2016, www.youtube.com/watch?v=vdNb9amv6dl.

C. Publications

Goeres, Darla M., et al. "Statistical assessment of a laboratory method for growing biofilms." *Microbiology* 151.3 (2005): 757-762.

Goeres, Darla M., et al. "Development, standardization, and validation of a biofilm efficacy test: The single tube method." *Journal of Microbiological Methods* 165 (2019): 105694.

Johnson, Erick, Theodore Petersen, and Darla M. Goeres. "Characterizing the Shearing Stresses within the CDC Biofilm Reactor Using Computational Fluid Dynamics." *Microorganisms* 9.8 (2021): 1709.

Buckingham-Meyer, Kelli, Darla M. Goeres, and Martin A. Hamilton. "Comparative evaluation of biofilm disinfectant efficacy tests." *Journal of microbiological methods* 70.2 (2007): 236-244.

Gomes, Inês B., et al. "Standardized reactors for the study of medical biofilms: a review of the principles and latest modifications." *Critical reviews in biotechnology* 38.5 (2018): 657-670.

Mendez, Ellen, et al. "The use of a CDC biofilm reactor to grow multi-strain *Listeria monocytogenes* biofilm." *Food Microbiology* 92 (2020): 103592.

Lindsay, D., et al. "The CDC biofilm bioreactor is a suitable method to grow biofilms, and test their sanitiser susceptibilities, in the dairy context." *International Dairy Journal* 126 (2022): 105264.

Williams, Dustin L., and Roy D. Bloebaum. "Observing the biofilm matrix of *Staphylococcus epidermidis* ATCC 35984 grown using the CDC biofilm reactor." *Microscopy and Microanalysis* 16.2 (2010): 143-152.

Ausbacher, Dominik, et al. "Paired methods to measure biofilm killing and removal: a case study with Penicillin G treatment of *Staphylococcus aureus* biofilm." *Letters in applied microbiology* 66.3 (2018): 231-237.